



4-H Embryology

How to open and pack up the 4-H Incubators


Open the incubator box from the top side that says, “Open This End”. The box is sealed with Velcro to keep the box lid closed.

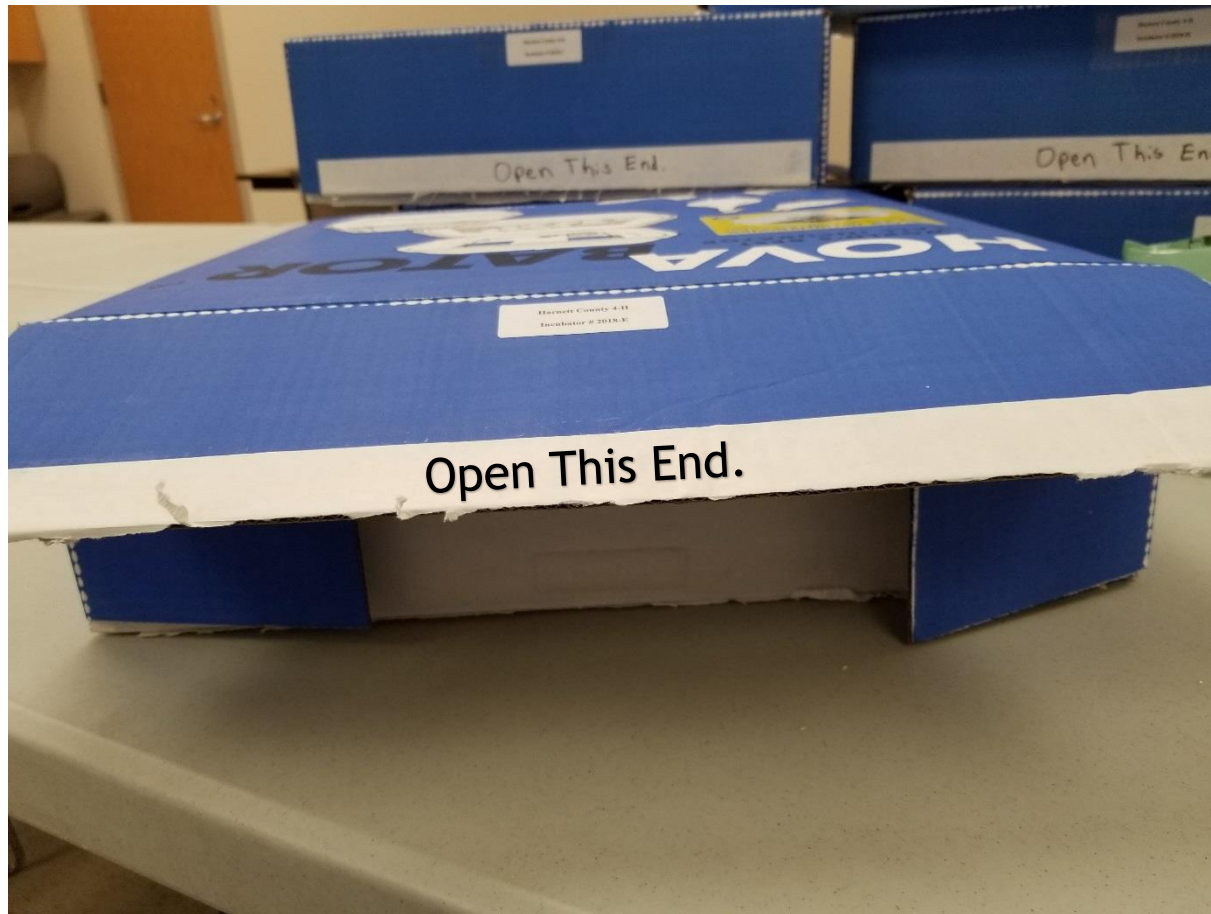


Pull incubator out from the sides.

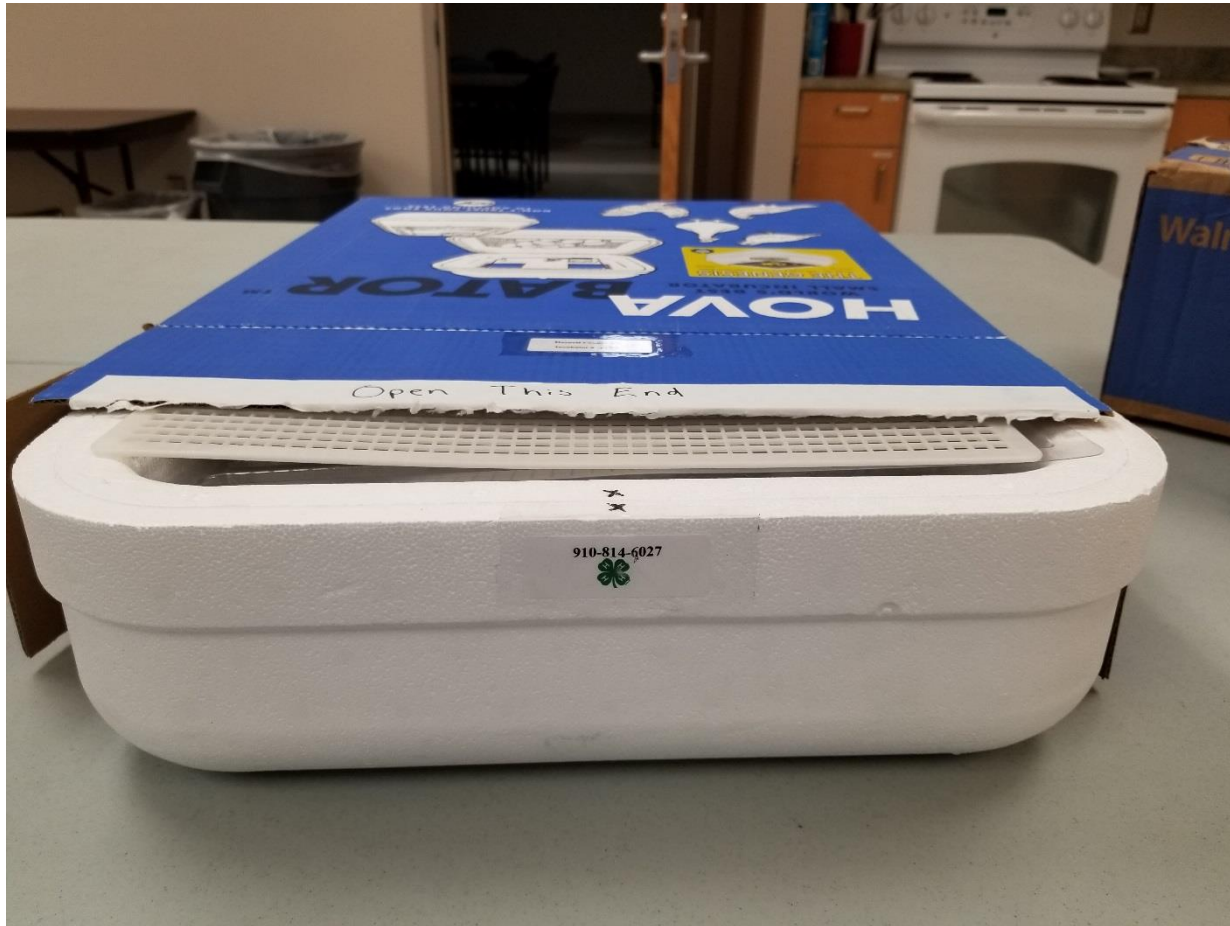

Please note that the plastic floor and clear plastic liner are on top of the incubator.

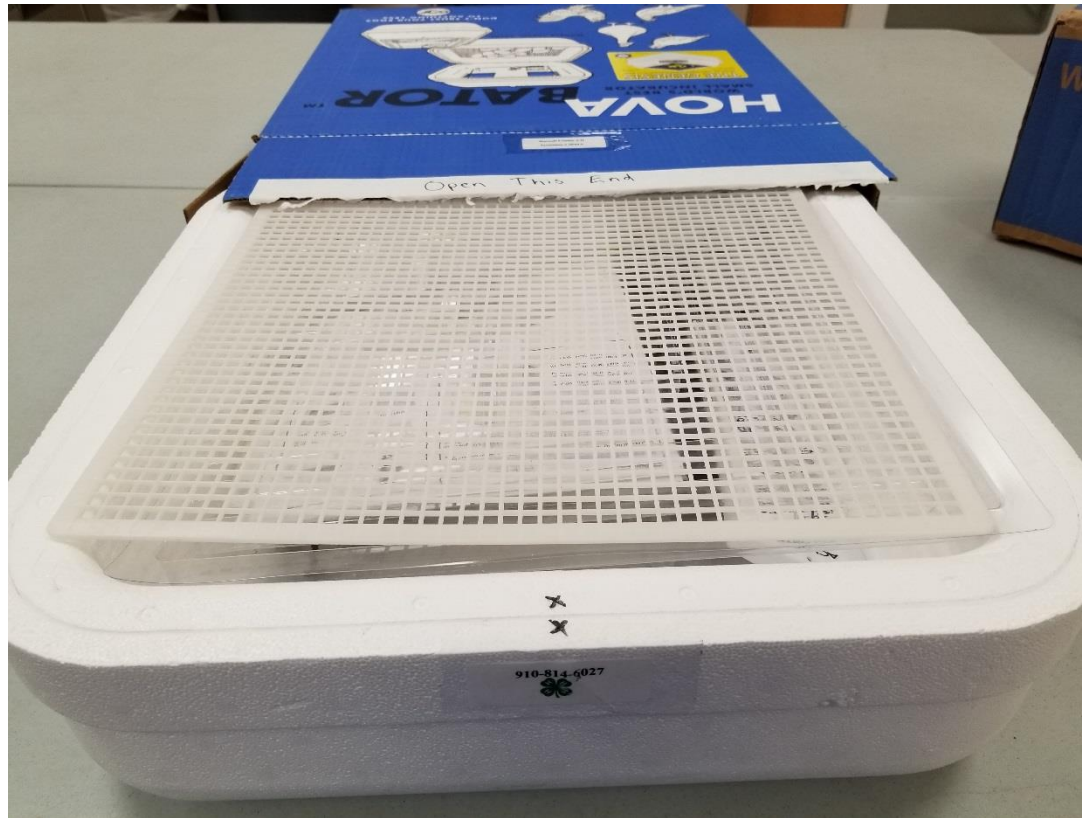


Remove the plastic floor and plastic liner and you will see thermometer, incubator information and hand washing poster.

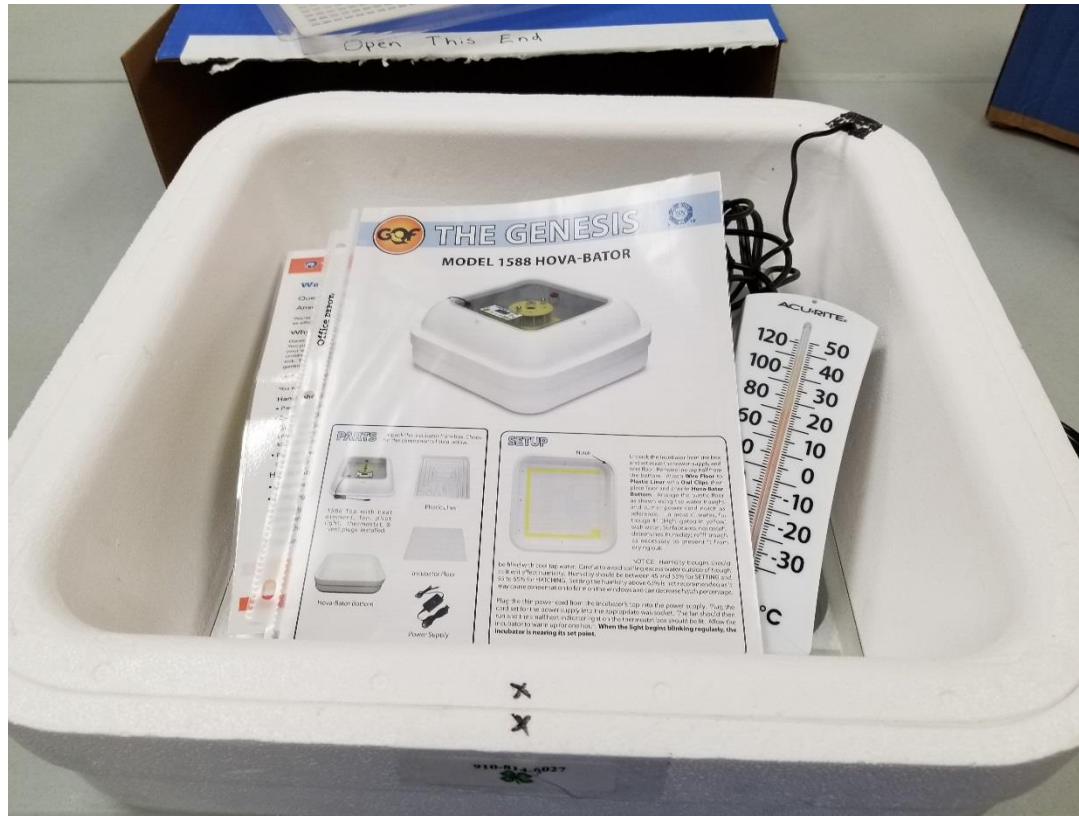

Once you remove the papers and thermometer, you have the electric cord.



Please notice that the wire comes through the hole in the back of the incubator. It has been marked in a black square.

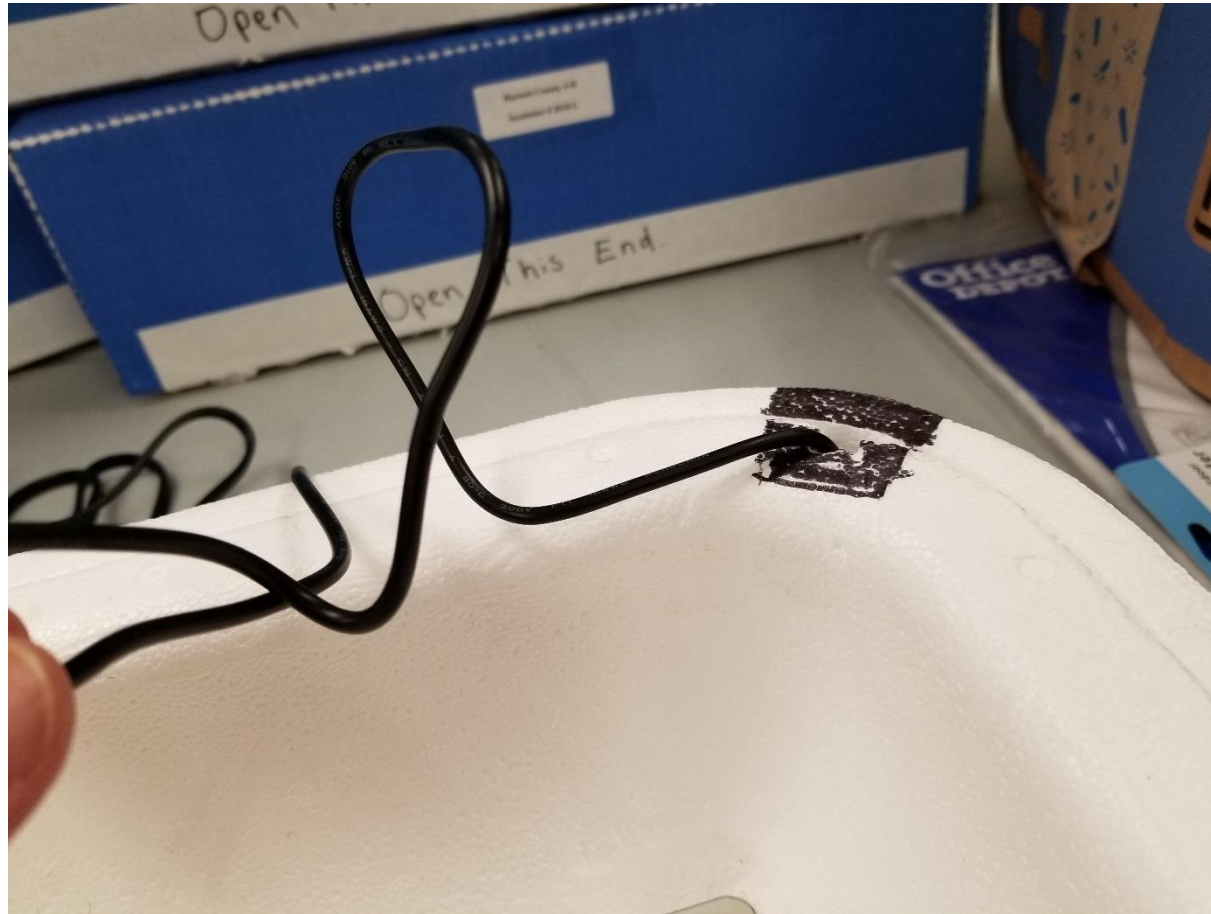

Remove the top of the incubator and place the clear plastic liner using the numbers to align it correctly. Add water to the liner in section 1 only till full.



Now add the plastic floor on top of the plastic liner.



Place the top on the incubator with the logos facing the front of your classroom for everyone to see and plug in.




After hatching, please wash out the bottom of the incubator, the plastic liner and plastic floor. The top can be wiped down if needed with a damp paper towel.




After everything is dry, pack top of incubator on bottom of incubator. Please make sure power cord come through hole in back.




Add thermometer and paperwork.

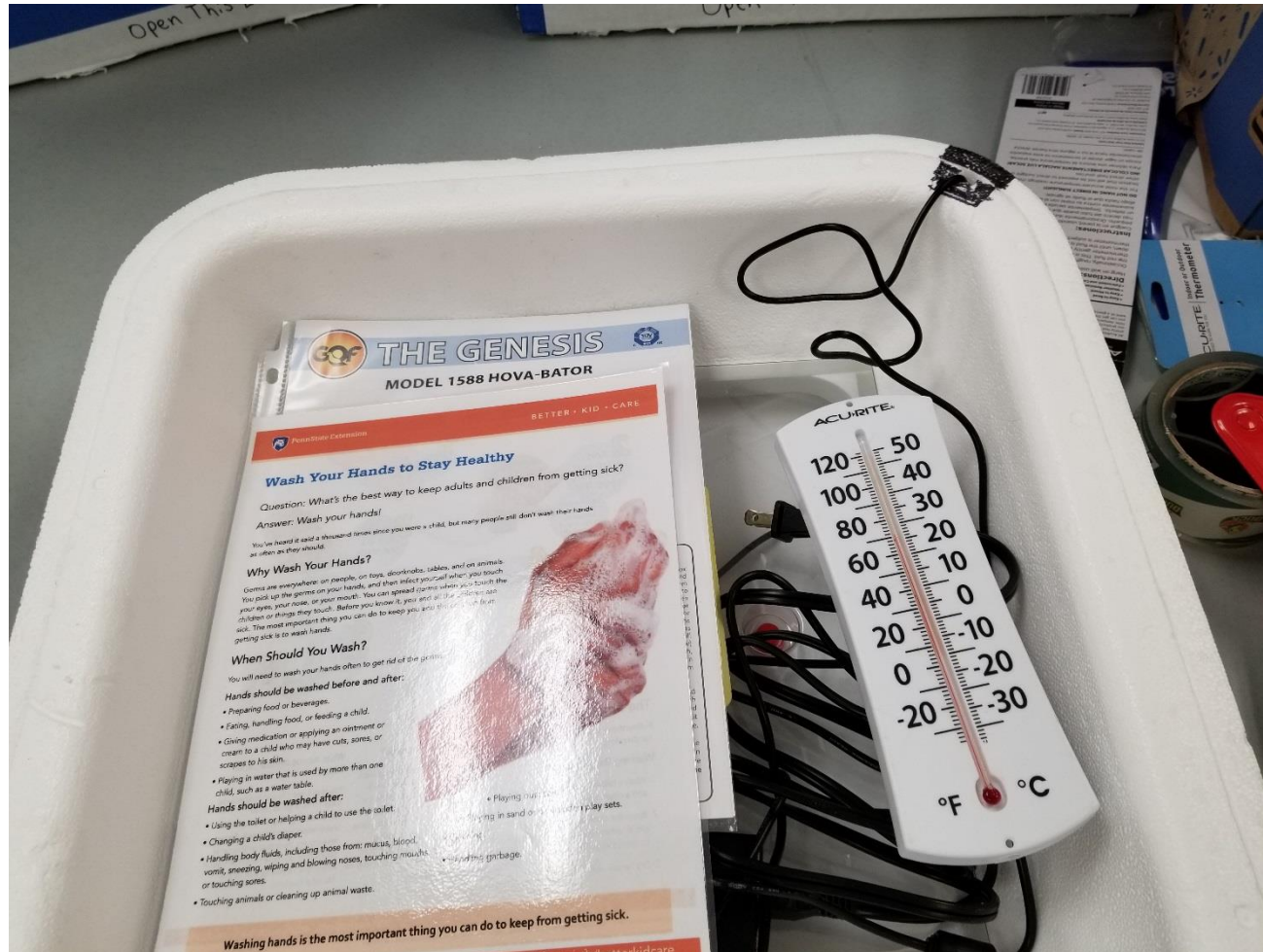

Place plastic liner and plastic floor on top of paperwork and insert to the box as shown below.

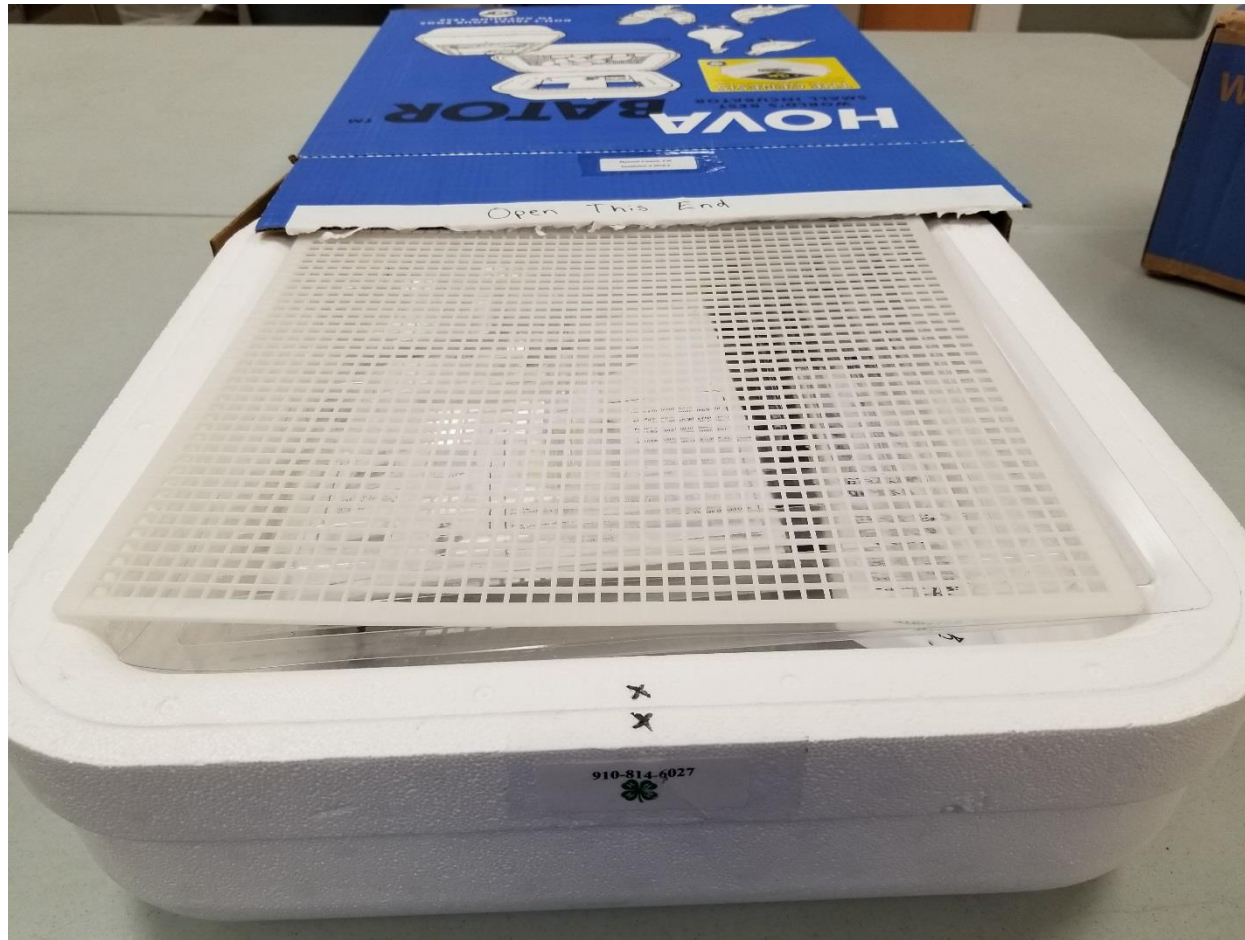

Push carefully into box.

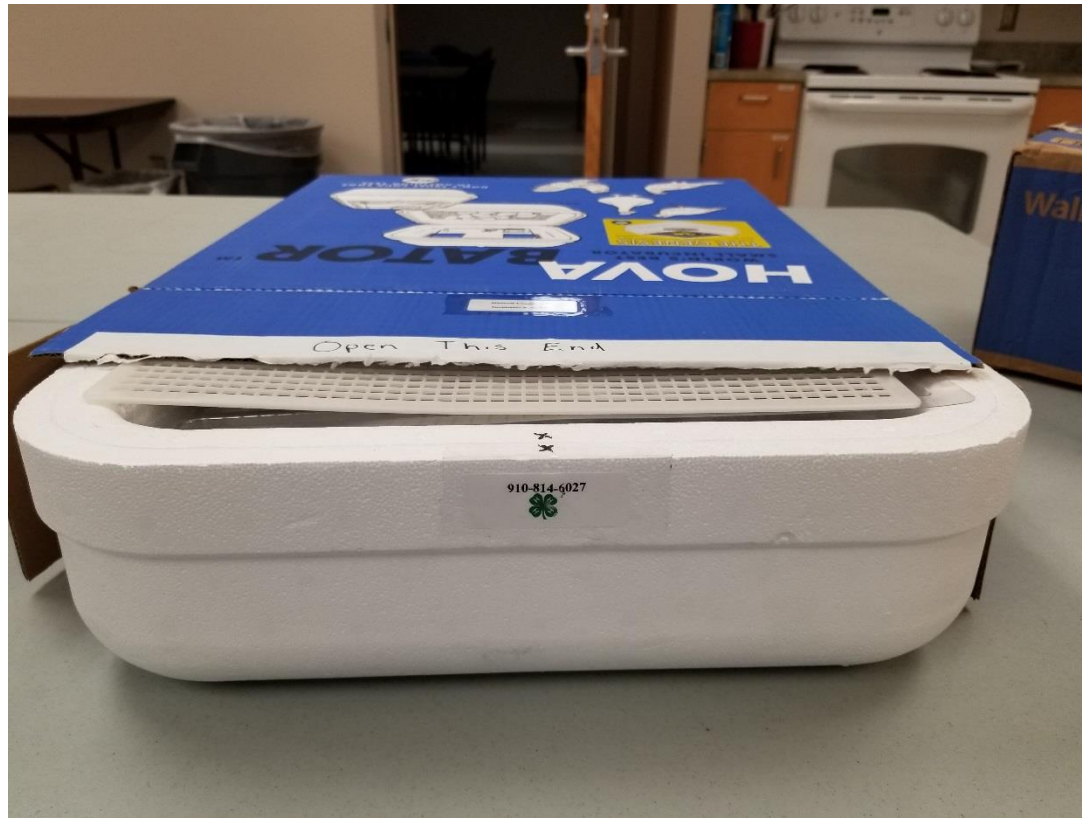

Fold white section on bottom first, then fold the two blue flaps. Then push top closed. Velcro will keep lid shut.

